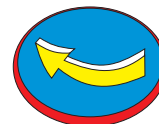
SSH

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Contents

Contents	3	Overview	
Warning and Safety Precautions	4	General View of the Scale	64
Getting Started	5	Description of the Keys	65
Operating Design	12	Specifications	66
Configuring the Scale		Dimensions (Scale Drawings)	68
Setting Parameters in the Operating Menu	17	Accessories (Options)	70
Scale Operating Menu (Overview)	20	CE Marking	72
Operating the Scale		Index	74
Basic Weighing Function	24		
Calibration/Adjustment	27		
Application Programs			
Counting	29		
Toggle between Weight Units	32		
Second Tare Memory	36		
Weighing in Percent	40		
Calculation by a Factor	43		
Totalizing	46		
Gross/Net Toggling	49		
Averaging	51		
Additional Functions	53		
Data Output Functions	55		
Printing a Data Record	55		
Error Codes	59		
Care and Maintenance	62		
Instructions for Recycling	63		

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This scale has been constructed in accordance with the European Directives as well as international regulations and standards for operation of electrical equipment, electromagnetic compatibility, and stipulated safety requirements.

Read these operating instructions thoroughly before using your scale to prevent damage to the equipment. Keep these instructions in a safe place.

Follow the instructions below to ensure safe and trouble-free operation of your scale:

- ⚠ Do not use this scale in a hazardous area/location
- ⚠ Make sure that the voltage rating printed on the AC adapter is identical to your local line voltage
- The scale housing is IP54-protected against dust deposits and water splashes
- The display and control unit is IP40-protected against penetration of solid foreign objects
- Connect only SCALTEC accessories and options, as these are optimally designed for use with your SSH scale
- Protect the AC adapter and the display and control unit from contact with liquids

When cleaning your scale, make sure that no liquids enters the scale housing; use only a slightly moistened cloth to clean the scale.

Do not open the scale housing. If the seal is broken, this will result in forfeiture of all claims under the manufacturer's warranty.

In case you have any problems with your scale:

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Warranty

Do not miss out on the benefits of our extended warranty. Please complete the warranty registration card and return it to the address listed below (see supplement).

In the unlikely event that your balance should require servicing within the 3-year warranty period, please contact your local SCALTEC dealer or:

SCALTEC INSTRUMENTS GmbH
Rudolf-Diesel-Strasse 1
37308 Heiligenstadt, Germany
Phone: (0)3606-6701-0
Fax: (0)3606-6701-60

Storage and Shipping Conditions

Do not expose the scale to shocks, vibrations, moisture or extreme temperatures.

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Unpacking the Scale

- After unpacking the scale, check it immediately for any visible damage as a result of rough handling during shipment.
- Note: The display and control unit is attached to the weighing platform via a cable.
- If you see any sign of damage, proceed as directed in the chapter entitled "Care and Maintenance," under the section on "Safety Inspection."

Save the box and all parts of the packaging until you have successfully installed your scale. Only the original packaging provides the best protection for shipment. Before packing your scale, unplug all connected cables to prevent damage.

Equipment Supplied

The equipment supplied includes the components listed below:

- Scale with attached display and control unit
- Load pan
- AC adapter
- Installation and operating instructions

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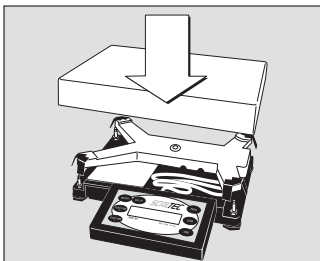
Installation Instructions

The SCALTEC scales are designed to provide reliable weighing results under normal ambient conditions. When choosing a location to set up your scale, observe the following so that you will be able to work with added speed and accuracy:


- Set up the scale on a stable, even surface
- Avoid placing the scale in close proximity to a heater or otherwise exposing the scale to heat or direct sunlight
- Protect the scale from drafts that come from open windows or doors
- Avoid exposing the scale to extreme vibrations during weighing
- Protect the scale from aggressive chemical vapors
- Do not expose the scale to extreme moisture over long periods

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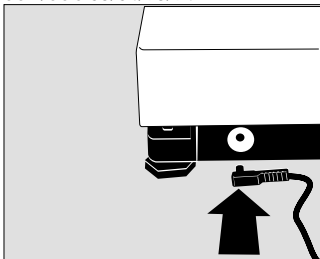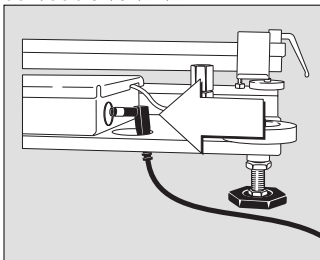
Connecting the Scale to AC Power

- Check the voltage rating and the plug design

Use only:

- Original SCALTEC AC adapters
- AC adapters with a registered approval rating from a national testing laboratory
- See the chapter entitled "Accessories" for information on using an external rechargeable battery pack with your scale
- Insert the right-angle plug into the jack and then tighten the screws
- Then insert the plug of the AC adapter into a wall outlet (mains)

SSH/SSC 81/82/91/92:


Information on Radio Frequency Interference

Warning!

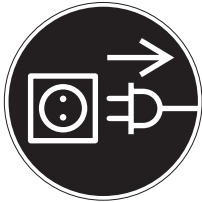This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause interference to radio communications. It has been tested and found to comply with the limits for a Class A computing device pursuant to Subpart J of Part 15 of FCC rules, which are designed to provide reasonable protection against such interference, when operated in a commercial environment. Operation of this equipment in a residential area is likely to cause interference, in which case the user, at his own expense, will be required to take whatever measures may be required to correct the interference.

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
Connecting Electronic Peripheral Devices

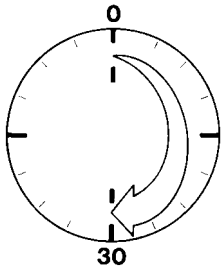- Make absolutely sure to unplug the scale from AC power before you connect or disconnect a peripheral device (printer or PC) to or from the interface port.

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
Warmup Time

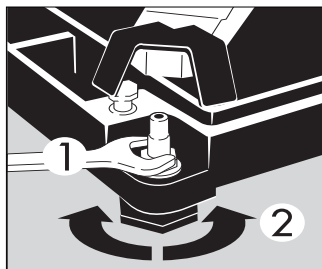To deliver exact results, the scale must warm up for at least 30 minutes after initial connection to AC power. Only after this time will the scale have reached the required operating temperature.

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
Leveling the Scale

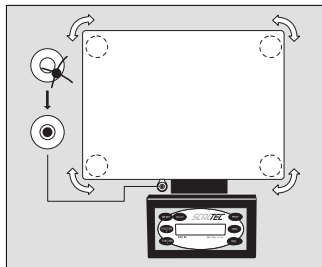- Remove the load plate
- Loosen the lock nuts on all leveling feet (use a 13 mm open-end wrench if necessary)
- Turn the leveling feet until the scale is level
- Avoid unintentional changes in the adjustment: Tighten the lock nuts with an open-end wrench and then replace the load plate

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Leveling the Scale (for SSH9. models only)

- Using the 4 foot screws, level the scale so that the air bubble is centered within the level indicator

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The scale consists of a weighing cell and a display and control unit. In addition to the choice of power supply (via AC adapter or external rechargeable battery pack), your scale also has an interface port for connecting peripheral devices, such as a printer, computer, universal remote control switch, etc.

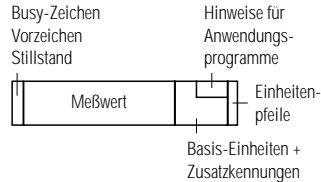
The display and control unit is affixed to the weighing platform. Operation of the scale is simple and uniform.

Keys

The scale can be operated using either the keys on the display and control unit or a connected computer (PC). Some of the keys activate different functions, depending on how long they are pressed ("press briefly" = < 2 sec.; "press and hold" = ≥ 2 sec.).

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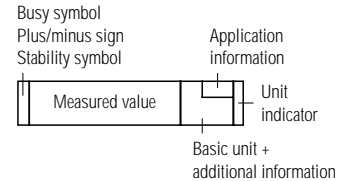
Display

The display is divided into 5 sections:

Busy Symbol, Plus/Minus Sign, Stability Symbol

If the symbol displayed here is

- a triangle ▲, this indicates that the scale is performing a function (busy symbol)
- a plus or minus sign (+ or –), this applies to the value displayed (this also applies to calculated values; e.g., during counting)
- a flashing triangle ▼, this indicates that the calculated values are not valid for legal metrology

Measured Value

When the scale is in the weighing mode, the weight, calculated values and application parameters are displayed here. In the setup mode, the menu code numbers are shown here.

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Basic Unit and Additional Information

When the scale has stabilized, the weight unit is displayed here (g or kg). The calculated values (e.g., % or pcs) are also displayed.

Application Information

This section shows information concerning the current application program; e.g., the reference quantity for counting.

Input Functions

When the scale is in the setup mode, you can select parameters from lists. The lists are contained in a menu, which is divided into 3 levels.

To set parameters: Turn the scale off and back on again. While all segments are displayed, press [TARE] briefly.

To navigate within a menu level: Press [TARE] briefly; when you reach the last menu item, the first is shown again.

To change menu levels: Press [PRINT] briefly.

To confirm the selected parameters on all 3 levels: Press and hold the [PRINT] key.

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To store parameter settings and exit the menu:

Press and hold the [TARE] key

Parameter settings are stored in the non-volatile memory. When you turn on the scale, the last parameters used are active.

To exit the menu without storing any new settings:

Press [ON/OFF]

Output Functions

Your scale is equipped with a data interface for connecting your choice of the following:

- Printer
- PC
- Universal remote control switch

Printer

You can configure a variety of options for data output to a printer.

You can have a printout generated automatically or only when the [PRINT] key is pressed. You can make the print function dependent on or independent of stability parameters, and define whether data ID codes will be included in the printout or not.

The print functions are described in detail under "Data Output Functions" in the chapter entitled "Operating the Scale."

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Data Interface

Instead of a printer, you can connect a different peripheral device, such as a computer, to the data interface.

You can use a connected computer to control the scale. Communication between the scale's data interface and a PC is in the form of request and response messages. Not all request telegrams will activate response messages.

See "Data Output Functions" in the chapter entitled "Operating the Scale" for a detailed description.

Error Messages

Error messages are displayed for 2 seconds. The format is as follows:

- Sequence error: "E" + 2 digits
- Hardware error: "E" + 3 digits

The section entitled "Error Messages" contains detailed information.

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Setting the Parameters (Menu)

Purpose

You can configure your scale to meet individual requirements by selecting from the parameters available in the menu.

Features

The parameters are divided into the following groups (1st menu level):

- 1 Scale functions
- 2 Application programs
- 3 Application parameters
- 5 Data interface
- 6 Print for weighing
- 7 Print for application
- 8 Extra functions
- 9 Reset menu to factory-set parameters

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Factory-Set Parameters

The factory-set parameters are marked by an "o" in the list starting on page 20.

Preparation

- Turn off the scale: Press [ON/OFF]
- Turn on the scale: Press [ON/OFF]; while all segments are displayed, press [TARE] briefly
- > Measured value line: ! (1st menu level)
- To navigate within a menu level: Press [TARE]; when you reach the last menu item, the first is shown again
- To select the next sublevel within a group: Press [PRINT] briefly
- To return to the higher menu level: Press [PRINT]
- To confirm selected parameters: Press and hold [PRINT] for more than two seconds
- > "o" indicates the selected menu code
- To store parameter settings and exit the menu: Press and hold the [TARE] key
- To exit the menu without storing any new settings: Press [ON/OFF]
- > Restart the application

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
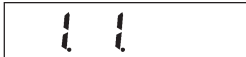
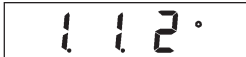
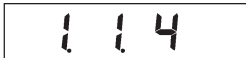
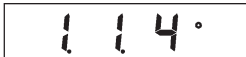
Example

Adapt the scale to unstable ambient conditions: Menu code 1.1.4

Step	Key	Display
1. Turn the scale off	[ON/OFF]	
2. Turn the scale back on; while all segments are displayed	[ON/OFF] [TARE] briefly	 
3. Confirm 1st menu level (scale functions)	[PRINT]	
4. Confirm "Adapt filter" group (2nd menu level)	[PRINT]	
5. Menu level 3: Select the desired menu item	[TARE] repeatedly	
6. Confirm selected item	[PRINT] for 2 seconds	
7. Set other parameters, if desired	[PRINT], [TARE]	
8. Saves changes and exit menu	[TARE] for 2 seconds	

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


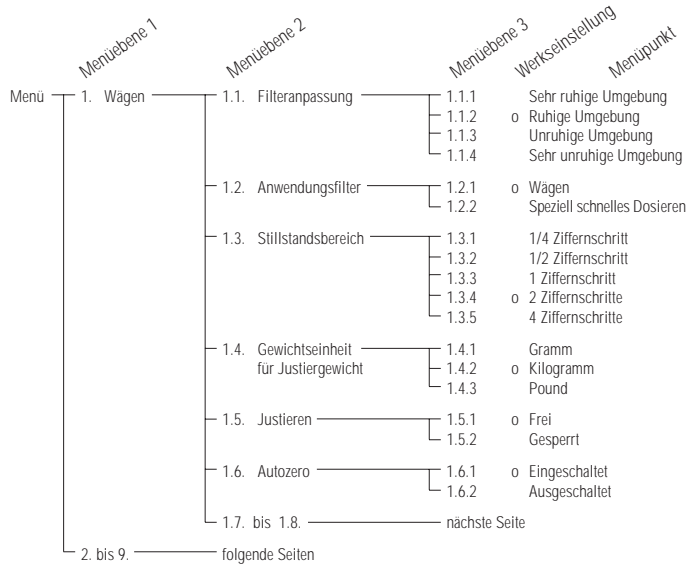
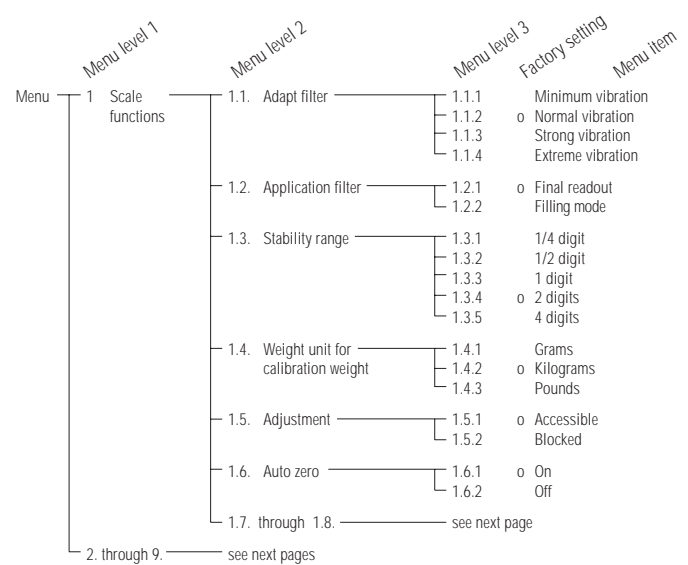
	Menu level 1	Menu level 2	Menu level 3	Factory setting	Menu item
Menu	1. Scale functions	1.1. through 1.6.	see previous page		
		1.7. Weight unit 1	1.7.2	Grams	
			1.7.3	o Kilograms	
			1.7.4	Carats	
			1.7.5	Pounds	
			1.7.6	Ounces	
			1.7.7	Troy ounces	
			1.7.8	Hong Kong taels	
			1.7.9	Singapore taels	
			1.7.10	Taiwanese taels	
			1.7.11	Grains	
			1.7.12	Pennyweights	
			1.7.13	Milligrams	
			1.7.14	Parts per pound	
			1.7.15	Chinese taels	
			1.7.16	Mommes	
			1.7.17	Austrian carats	
			1.7.18	Tola	
			1.7.19	Baht	
			1.7.20	Mesghal	
			1.7.21	Tons	
			1.7.22	Pounds+ounces (non-decimal display)	
		1.8. Display accuracy 1	1.8.1	o Standard resolution	
			1.8.3	2x higher resolution (PolyRange)	
	2. Application programs	2.1. Program selection	2.1.1	Weighing	
			2.1.2	Toggleing weight units	
			2.1.3	Net-total (tare memory)	
			2.1.4	o Counting	
			2.1.5	Weighing in percent	
			2.1.7	Calculation by a factor	
			2.1.8	Totalizing	
			2.1.9	Gross/net toggleing	
			2.1.10	Tare memory with autom. container taring	
			2.1.11	Weighing in percent with difference display	
			2.1.12	Averaging	
	3. through 9.	see following pages			

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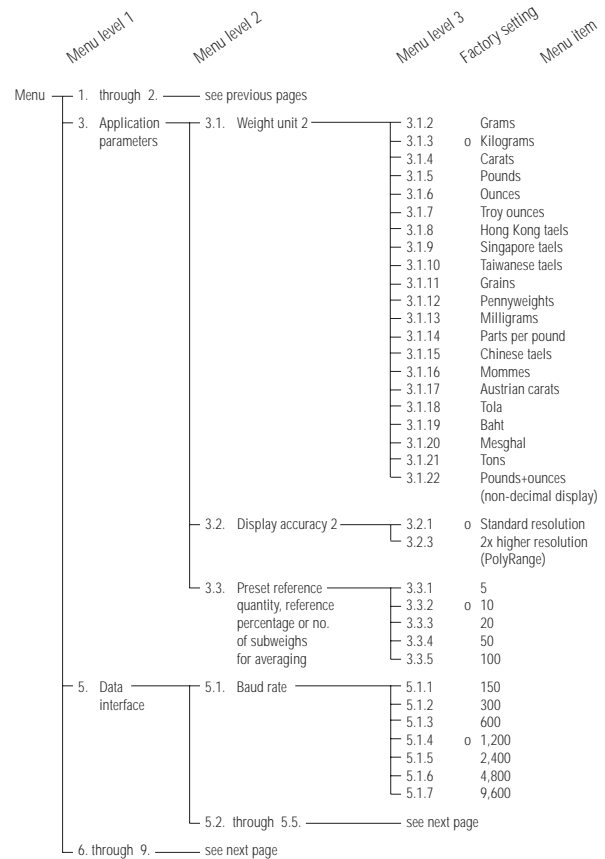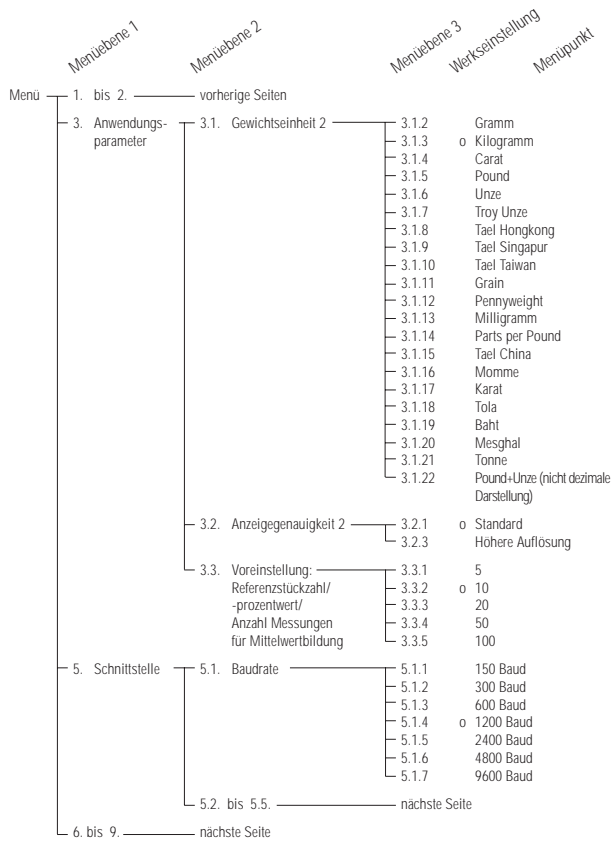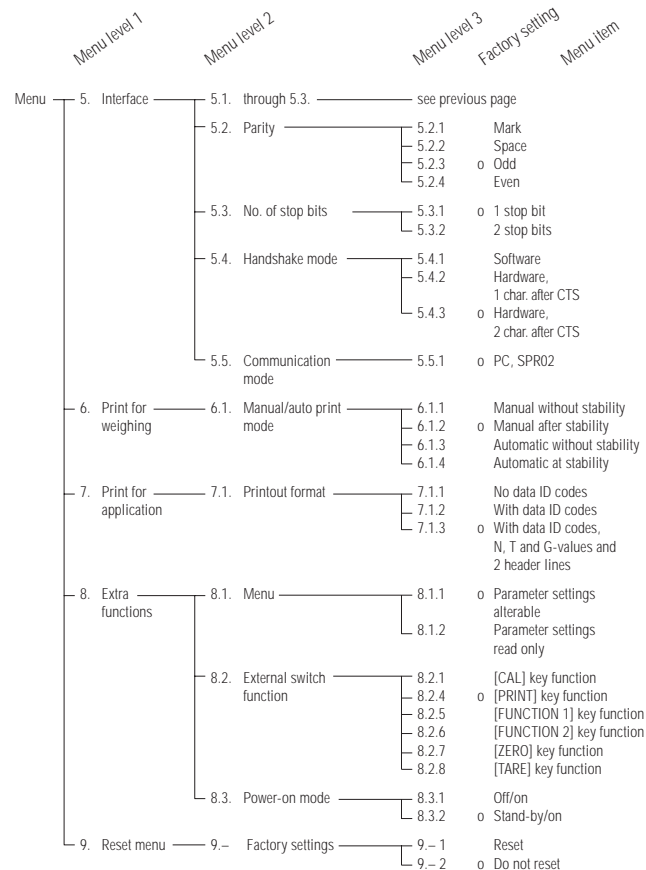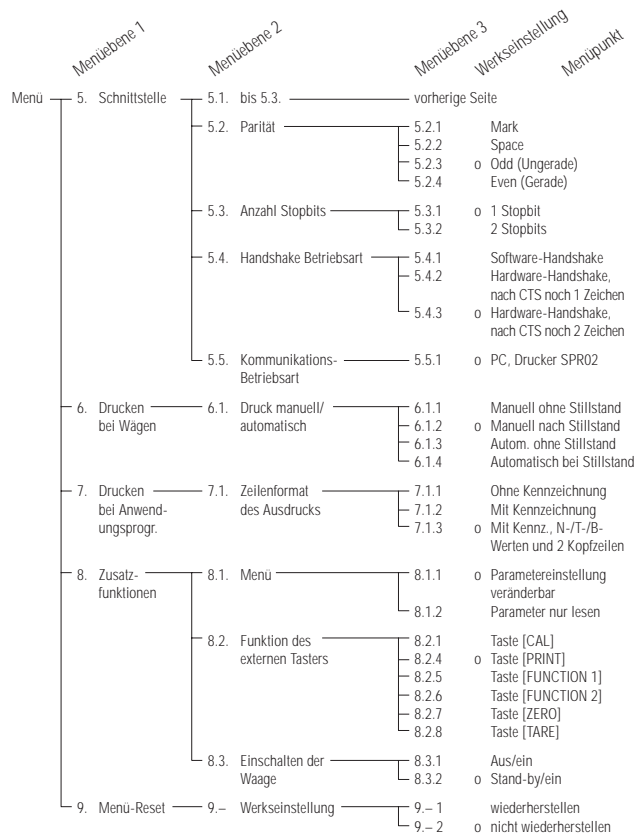Basic Weighing Function

Purpose

The basic weighing function can be used alone or in combination with an application program (Toggle between Weight Units, Counting, Weighing in Percent, etc.).

Features

- Zeroing the scale
Depending on prevailing ambient conditions, the display may not show a zero readout even though there is no load on the scale. If the weight shown is less than 2% of the scale's maximum weighing capacity, you can zero the scale.
- Taring the scale
Tare the scale with an empty container on the weighing pan to obtain a readout of the net weight after the container is filled.
- Printing weights

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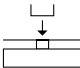
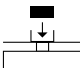
Preparation

- Turn on the scale: Press [ON/OFF]
- > All segments light up briefly
- To change configurations: See the chapter entitled "Configuring the Scale"
- To load factory-set configurations: See "Configuring the Scale," parameter 9 - 1

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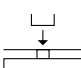
* The SCALTEC Service Center or your SCALTEC dealer can configure the scale to include 2 customer-specific lines on the printout. Software is available for setting this configuration with a PC.

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Calibration/Adjustment

Purpose

Calibration is the determination of the difference between the weight readout and the true weight (mass) of a sample. Calibration does not entail making any changes within the scale.

Adjustment is the correction of this difference between the measured value displayed and the true weight (mass) of the sample, or the reduction of the difference to an allowable level within the maximum permissible error limits.

Features

External adjustment can only be performed when

- there is no load on the scale,
- the scale is set to zero, and
- the internal signal is stable.

If these conditions are not met, an error message is displayed.

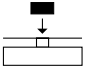
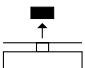
The weight on the scale must not differ from the nominal weight by more than 2%.

You can use g, kg or lb as the weight unit for calibration (! 4)

You can block external calibration of the scale (! 5)

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

Example

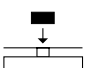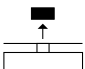Calibrate the scale

Step	Key (or instruction)	Display/Output
1. Turn on the scale	[ON/OFF]	0.000 kg
2. If necessary, zero the scale	[ZERO]	0.000 kg
3. Begin calibration Calibration weight is displayed without wt. unit	[CAL/CF]	+ 5.000
4. Place the indicated calibration weight on the scale (here: 5,000 g) After calibration, the calibration weight is displayed with wt. unit		5.000 + 5.000 kg
5. Remove the calibration weight		0.000 kg

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Counting

Purpose

With the Counting program you can determine the number of parts that have approximately the same average piece weight.

Available Features

- The reference sample quantity can be changed either in the operating menu or during weighing
- Store the current weight value to have it loaded as the preset reference sample quantity the next time you initialize the Counting application
- Re-initialize without quitting the program
- The average piece weight is automatically output via the data interface port after initialization, if the menu code for "Printout with data ID codes" is set
- Press [FUNCTION 2] to toggle the display between piece count and weight

Preparation

Configure the Counting application in the operating menu:

- If the scale is on, turn it off: Press [ON/OFF]
- Turn on the scale: Press [ON/OFF]; while all segments are lit, press [TARE]

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- Confirm the selection of Counting: Press [PRINT] for at least 2 sec.

> o indicates that this is the current menu code setting

- Set the next parameters: Press [PRINT]

- Select and confirm:

- Reference sample quantity:

3 3 1 5 pcs
3 3 2 10 pcs
3 3 3 20 pcs
3 3 4 50 pcs
3 3 5 100 pcs

see also the “Scale Operating Menu (Overview)” in the chapter entitled “Configuring the Scale”

- Save settings and exit the Setup menu: Press [TARE] for at least 2 sec.

Additional Functions

In addition to the basic functions (power off, zeroing, taring and printing), you can also access the following functions from this application:

- Initialize application ([FUNCTION 1] key)
- Delete initialization values ([CAL/CF] key; press for at least 2 sec.)
- Change reference sample quantity after deleting initialization value ([FUNCTION 2] key, press for at least 2 sec.)
- Display weight ([FUNCTION 2] key)
- Calibrate/adjust scale ([CAL/CF] key)

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Example

Determine an unknown piece count; store the weight on the scale as a reference sample quantity

Settings (changes in the factory settings required for this example):

Menu: Application program: Counting (2 / 4)

Step	Key (or instruction)	Display/Output
1. Turn on the scale	[ON/OFF]	
2. If necessary, zero the scale	[ZERO]	0.000 kg ^{kg}
3. Select the reference sample quantity (here: 50 pcs)	[FUNCTION 2] [FUNCTION 2]	rEF 20 (briefly) rEF 50 (briefly)
4. Place the reference sample quantity (50 pcs) on the scale (here: 0.930 kg)		+ 0.930 kg ^{kg}
5. Start the application; if the print format is set to include data ID codes, the following is printed	[FUNCTION 1]	+ 50 pcs ^{pcs}
6. Display weight	[FUNCTION 2]	wRef + 0.186 kg
7. Display quantity	[FUNCTION 2]	+ 0.930 kg ^{kg}
8. Weigh uncounted parts (here: 174 pcs)		+ 50 pcs ^{pcs}
9. Print total piece count	[PRINT]	+ 174 pcs ^{pcs}
10. Unload the scale		EISENMUELLER KASSEL Qnt + 174 pcs
11. Repeat the procedure starting from Step 8, if desired.		0 pcs ^{pcs}

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Toggle between Weight Units

Purpose

With this application program you can toggle the display of a weight value back and forth between two weight units.

Available Features

- Toggling the displayed weight
- Other features as for the basic weighing function

Preparation

Configure the Toggle weight units application in the operating menu:

- If the scale is on, turn it off: Press [ON/OFF]
- Turn on the scale: Press [ON/OFF]; while all segments are lit, press [TARE]
- Select the "Toggle weight units" menu item: Press [TARE], [PRINT], and [PRINT]; press [TARE] repeatedly, if necessary
2 1 2 Toggle weight units
- Confirm selection: Press [PRINT] for more than 2 sec.
> o indicates that this is the current menu code setting
- Set the next parameters: Press [PRINT]

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● Select and confirm:

- Weight unit 1: see next page (1 7 x)
- Display resolution 1:
 - ! B ! Standard resolution
 - ! B 3 2x higher resolution (PolyRange)
- Weight unit 2: see next page (3 1 x)
- Display resolution 2:
 - 3 2 ! Standard resolution
 - 3 2 3 2x higher resolution (PolyRange)

see also the "Scale Operating Menu (Overview)" in the chapter entitled "Configuring the Scale"

● Save settings and exit the Setup menu: Press [TARE] for at least 2 sec.

Additional Functions

In addition to the functions for:

- power off ([ON/OFF] key),
- zeroing ([ZERO] key),
- taring ([TARE] key), and
- printing ([PRINT] key),

you can also access the following functions from this application:

- toggle weight unit ([FUNCTION 2] or [FUNCTION 1] key),
- calibration/adjustment ([CAL/CF] key)

Menüpunkt	Einheit	Umrechnung	Druck
(7 2) (3 2)	Gramm	1,0000000000	g
(7 3) (3 3)	Kilogramm	0,0010000000	kg
(7 4) (3 4)	Carat	5,0000000000	ct
(7 5) (3 5)	Pound	0,00220462260	lb
(7 6) (3 6)	Ounce	0,03527396200	oz
(7 7) (3 7)	Troy ounce	0,03215074700	ozt
(7 8) (3 8)	Tael Hongkong	0,02671725000	tlh
(7 9) (3 9)	Tael Singapur	0,02645544638	tls
(7 10) (3 10)	Tael Taiwan	0,02666666000	tlt
(7 11) (3 11)	Grain	15,43235835000	GN
(7 12) (3 12)	Pennyweight	0,64301493100	dwt
(7 13) (3 13)	Milligramm	1000,0000000000	mg
(7 14) (3 14)	Parts per Pound	1,12876677120	/lb
(7 15) (3 15)	Tael China	0,02645547175	tlc
(7 16) (3 16)	Momme	0,26670000000	mom
(7 17) (3 17)	Karat	5,0000000000	K
(7 18) (3 18)	Tola	0,08573333810	tol
(7 19) (3 19)	Baht	0,06578947437	bat
(7 20) (3 20)	Mesghal	0,21700000000	MS
(7 21) (3 21)	Tonne	0,00000100000	T
(7 22) (3 22)	lb/oz	0,03527396200	o

Menu code	Unit	Conversion factor	Printout
(7 2) (3 2)	Grams	1.00000000000	g
(7 3) (3 3)	Kilograms	0.00100000000	kg
(7 4) (3 4)	Carats	5.00000000000	ct
(7 5) (3 5)	Pounds	0.00220462260	lb
(7 6) (3 6)	Ounces	0.03527396200	oz
(7 7) (3 7)	Troy ounces	0.03215074700	ozt
(7 8) (3 8)	Hong Kong taels	0.02671725000	tlh
(7 9) (3 9)	Singapore taels	0.02645544638	tls
(7 10) (3 10)	Taiwanese taels	0.02666666000	tlt
(7 11) (3 11)	Grains	15.43235835000	GN
(7 12) (3 12)	Pennyweights	0.64301493100	dwt
(7 13) (3 13)	Milligrams	1000.0000000000	mg
(7 14) (3 14)	Parts per pound	1.12876677120	/lb
(7 15) (3 15)	Chinese taels	0.02645547175	tlc
(7 16) (3 16)	Mommes	0.26670000000	mom
(7 17) (3 17)	Austrian carats	5.00000000000	K
(7 18) (3 18)	Tola	0.08573333810	tol
(7 19) (3 19)	Baht	0.06578947437	bat
(7 20) (3 20)	Mesghal	0.21700000000	MS
(7 21) (3 21)	Tons	0.00000100000	T
(7 22) (3 22)	lb/oz	0.03527396200	o

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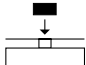
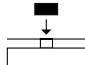
Example

Toggle the display from kilograms [kg] (1st unit) to pounds [lb] (2nd unit)

Settings (changes in the factory settings required for this example):

Menu: Application program: Toggle between weight units (2 | 2)

Menu: Weight unit 2: pounds (3 | 5)

Step	Key (or instruction)	Display/Output
1. Turn on the scale	[ON/OFF]	
2. Place weight on scale (here: 2.295 kg)		
3. Toggle to weight unit 2: Pounds [lb]	[FUNCTION 2]	
4. Print weight	[PRINT]	EISENMUELLER KASSEL G + 5.060 lb
5. Toggle to weight unit 1: kilograms [kg]	[FUNCTION 2]	

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Second Tare Memory

Purpose

With this application program you can store the weight on the scale as a tare compensation weight.

Available Features

- Press [FUNCTION 1] to store the weight on the scale in the second tare memory
- NET is displayed next to the net weight readout if the scale has been tared or if a tare value is stored in the second tare memory (**N** in the printout indicates that the scale was tared by pressing the [TARE] key; **N 1** indicates that there is data in the second tare memory)
- You can have a series of containers tared in sequence, if the scale is unloaded to under 30% of the previous weight after measurement and then loaded with 70% – 130% of the previous weight (i.e., the next container)
- Press [CAL/CF] for at least 2 seconds to delete the value from the second tare memory
- Press [FUNCTION 2] to toggle between weight unit 1 and weight unit 2

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Preparation

Configure the Second tare memory application in the operating menu:

- If the scale is on, turn it off: Press [ON/OFF]
- Turn on the scale: Press [ON/OFF]; while all segments are lit, press [TARE]
- Select the "Second tare memory" menu item: Press [TARE], [PRINT], and [PRINT]; press [TARE] repeatedly, if necessary
 - 2 1 3 Second tare memory
 - 2 1 10 Second tare memory with automatic container taring
- Confirm selection of Second tare memory: Press [PRINT] for at least 2 sec.
- > o indicates that this is the current menu code setting
- Set the next parameters: Press [PRINT]
- Select and confirm:
 - Weight unit 1: (see also "Toggle between Weight Units")
 - 1 7 2 Grams
 -
 - 1 7 2 2 lb / oz
 - Weight unit 2: (see also "Toggle between Weight Units")
 - 3 1 2 Grams
 -
 - 3 1 2 2 lb / oz

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see also the “Scale Operating Menu (Overview)” in the chapter entitled “Configuring the Scale”

- Save settings and exit the Setup menu:
Press [TARE] for at least 2 sec.

Additional Functions

In addition to the basic functions (power off, zeroing, taring and printing), you can also access the following functions from this application:

- Store weight in second tare memory ([FUNCTION 1] key)
- Delete tare memory ([CAL/CF] key; press for at least 2 sec.)
- Toggle weight unit ([FUNCTION 2] key)

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Example

Fill containers with nearly identical weights ($\pm 30\%$)

Settings (changes in the factory settings required for this example): Menu:

Application program: Second tare memory with automatic container taring ($\text{2} \text{ : } \text{!} \text{ : } \text{!} \text{ : } \text{!}$)

Menu: Weight unit 2: Grams ($\text{3} \text{ : } \text{2}$)

Step	Key (or instruction)	Display/Output
1. Turn on the scale	[ON/OFF]	
2. If necessary, zero the scale	[ZERO]	0.000 kg
3. Place an empty container on the scale (here: 65 g)		+ 0.065 kg
4. Store the container weight in the second tare memory. If the print format is set to include data ID codes, the following is printed	[FUNCTION 1]	0.000 kg _{NET}
5. Fill the container (here: to 500 g)		N1 + 0.065 kg + 0.500 kg _{NET}
6. Display weight unit 2	[FUNCTION 2]	+ 500 g _{NET}
7. Display weight unit 1	[FUNCTION 2]	+ 0.500 kg _{NET}
8. Remove the filled container from the scale		- 0.065 kg _{NET}
9. Place the next empty container on the scale (here: 75 g)		0.000 kg _{NET} (automatic container taring is performed)
10. Repeat the procedure starting from Step 5.		

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Weighing in Percent

Purpose

This application program allows you to obtain weight readouts in percent which are in proportion to a reference weight.

Available Features

- Reference percentage can be changed either in the operating menu or during weighing
- Store the current weight value to have it loaded as the preset reference percentage the next time you initialize the Weighing in Percent application
- Optional auto zero function
- The reference sample weight is automatically output via the data interface port after initialization, if the menu code for "Printout with data ID codes" is set
- Press [FUNCTION 2] to toggle the display between percentage and weight

Preparation

Configure the Weighing in Percent application in the operating menu:

- If the scale is on, turn it off: Press [ON/OFF]
- Turn on the scale: Press [ON/OFF]; while all segments are lit, press [TARE]

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- Select the “Weighing in Percent” menu item:
Press [TARE], [PRINT], and [PRINT]; press [TARE] repeatedly, if necessary
2 1 5 Weighing in percent – without taring
2 1 1 1 Weighing in percent – with taring

- Confirm selection of Weighing in Percent: Press [PRINT] for at least 2 sec.

> o indicates that this is the current menu code setting

- Set the next parameters: Press [PRINT]

- Select and confirm:
 - Reference percentage at power-on:
 - 3 3 1 5 %
 - 3 3 2 10 %
 - 3 3 3 20 %
 - 3 3 4 50 %
 - 3 3 5 100 %

see also the “Scale Operating Menu (Overview)” in the chapter entitled “Configuring the Scale”

- Save settings and exit the Setup menu: Press [TARE] for at least 2 sec.

Additional Functions

In addition to the basic functions (power off, zeroing, taring and printing), you can also access the following functions from this application:

- Initialize application ([FUNCTION 1] key)
- Delete initialization values ([CAL/CF] key; press for at least 2 sec.)
- Change reference percentage after deleting initialization value ([FUNCTION 2] key; press for at least 2 sec.)
- Display weight ([FUNCTION 2] key)
- Calibrate/adjust scale ([CAL/CF] key)

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Example

Determine an unknown percentage; store the weight on the scale as a reference percentage

Settings (changes in the factory settings required for this example):

Menu: Application program: Weighing in percent (2 : 5)

Menu: Reference percentage 100 % (3 3 5)

Step	Key (or instruction)	Display/Output
1. Turn on the scale	[ON/OFF]	
2. If necessary, zero the scale	[ZERO]	0.000 kg
3. Place the reference weight on the scale (here: 2.295 kg = 100%)		+ 2.295 kg
4. Start the application; if the print format is set to include data ID codes, the following is printed	[FUNCTION 1]	+ 100.00 % Wxx% + 2.295 kg
5. Display weight	[FUNCTION 2]	+ 2.295 kg
6. Display percentage	[FUNCTION 2]	+ 100.00 %
7. Display the reference percentage	[FUNCTION 1] (at least 2 sec.)	rEF 100
8. Place an unknown weight on the scale (here: 3.225 kg)		+ 140.41 %
9. Print percentage	[PRINT]	Prc + 140.41 %
10. Unload the scale		0.00 %
11. Repeat the procedure starting with Step 8, if desired.		

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Calculation by a Factor

Purpose

With this application program you can calculate the weight of a sample by a given factor.

Available Features

- Calculate a weight value using one of the following factors: 0.25; 0.50; 0.75; 1.0; 1.5; 2.0; 2.5; 3.0; 3.5; 4.0; 4.5; 5.0
- A flashing triangle in the display indicates a calculated value
- Press the [CAL/CF] key for at least 2 seconds to delete the calculation factor
- Press [FUNCTION 2] to toggle between weight unit 1 and weight unit 2

Preparation

Configure the Calculation application in the operating menu:

- If the scale is on, turn it off: Press [ON/OFF]
- Turn on the scale: Press [ON/OFF]; while all segments are lit, press [TARE]
- Select the "Calculation" menu item: Press [TARE], [PRINT], and [PRINT]; press [TARE] repeatedly, if necessary
2 1 7 Calculation by a factor

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- Confirm selection of Calculation: Press [PRINT] for at least 2 sec.

> o indicates that this is the current menu code setting

- Set the next parameters: Press [PRINT]

- Select and confirm:

- Weight unit 1: (see also "Toggle between Weight Units")

1 7 2 Grams

...

1 7 2 2 lb/oz

- Weight unit 2: (see also "Toggle between Weight Units")

3 1 2 Grams

...

3 1 2 2 lb/oz

see also the "Scale Operating Menu (Overview)" in the chapter entitled "Configuring the Scale"

- Save settings and exit the Setup menu: Press [TARE] for at least 2 sec.

Additional Functions

In addition to the basic functions (power off, zeroing, taring and printing), you can also access the following functions from this application:

- Initialize application ([FUNCTION 1] key)
- Change the calculation factor after the application is initialized ([FUNCTION 1] key)
- Delete initialization values ([CAL/CF] key; press for at least 2 sec.)
- Toggle weight unit ([FUNCTION 2] key)
- Calibration/adjustment ([CAL/CF] key)

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Example

You have a formula requiring several different components for a total amount of 1,000 g. You want to use this recipe to make a 500-g batch (factor: 0.5). The scale shows the amounts called for in the formula, although only half the amounts are placed on the scale. This saves you having to recalculate the amount of each component.

Settings (changes in the factory settings required for this example):

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Totalizing

Purpose

This application program acts as a cumulative memory function.

Available Features

- The current weight is stored as a net value when the scale stabilizes
- A value indicated as negative (–) in the display is stored as a positive (+) value in the totalizing memory
- The data stored in the totalizing memory is automatically output via the data interface port after initialization, if the menu code for "Printout with data ID codes" is set
- The number of values in the totalizing memory is displayed (transaction counter)
- Press [FUNCTION 1] for at least 2 sec. to display the contents of the totalizing memory
- Press [CAL/CF] to clear the totalizing memory. The total is displayed for 1.5 seconds and printed
- You cannot toggle between weight units when there is data in the totalizing memory

Preparation

Configure the Totalizing application in the operating menu:

- If the scale is on, turn it off: Press [ON/OFF]
- Turn on the scale: Press [ON/OFF]; while all segments are lit, press [TARE]

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- Select the "Totalizing" menu item:
Press [TARE], [PRINT], and [PRINT]; press [TARE] repeatedly if necessary
2 1 8 Totalizing
- Confirm selection of Totalizing: Press [PRINT] for at least 2 sec.
> o indicates that this is the current menu code setting
- Set the next parameters: Press [PRINT]
- Select and confirm:
 - Weight unit 1: (see also "Toggle between Weight Units")
1 7 2 Grams
 - ...
 - 1 7 2 2 lb/oz

see also the "Scale Operating Menu (Overview)" in the chapter entitled "Configuring the Scale"
- Save settings and exit the Setup menu: Press [TARE] for at least 2 sec.

Additional Functions

In addition to the basic functions (power off, zeroing, taring and printing), you can also access the following functions from this application:

- Start application ([FUNCTION 1] key)
- Display value in totalizing memory ([FUNCTION 1] key; press for at least 2 sec.)
- Clear totalizing memory ([CAL/CF] key; press for at least 2 sec.)

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Example

Totalize weight values

Settings (changes in the factory settings required for this example):

Menu: Application program: Totalizing (2 : 8)

Step	Key (or instruction)	Display/Output
1. Turn on the scale	[ON/OFF]	
2. If necessary, zero the scale	[ZERO]	0.000 kg
3. Place sample on the scale (here: 380 g)		+ 0.380 kg
4. Store value in memory; if the print format is set to include data ID codes, the following is printed	[FUNCTION 1]	+ 0.380 kg ¹ g + 0.380 kg
5. Place the next sample on the scale (here: 575 g)		+ 0.575 kg ¹
6. Store value in memory	[FUNCTION 1]	+ 0.575 kg ²
7. Display the total in memory (total is also printed)	[FUNCTION 1] for at least 2 sec.	±0.955 kg ² (briefly) sum + 0.955 kg
8. Clear totalizing memory and print total	[CAL/CF] at least 2 sec.	sum + 0.955 kg
9. Repeat the procedure starting from Step 4, if desired.		

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Gross/Net Toggling

Purpose

With this application program you can toggle between net and gross values.

Available Features

- Toggle the display between gross and net weights when there is data in the tare memory
- Press [FUNCTION 2] to toggle between weight unit 1 and weight unit 2

Preparation

Configure the Gross/net toggling application in the operating menu:

- If the scale is on, turn it off: Press [ON/OFF]
- Turn on the scale: Press [ON/OFF]; while all segments are lit, press [TARE] briefly
- Select the "Gross/net toggling" menu item: Press [TARE], [PRINT], and [PRINT]; press [TARE] repeatedly, if necessary
2 19 Gross/net toggling
- Confirm selection of Gross/net toggling: Press [PRINT] for at least 2 sec.
- > 0 indicates that this is the current menu code setting
- Set the next parameters: Press [PRINT]

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Averaging

Purpose

Use this program to determine the weights of unstable samples (e.g., live animals) or to determine weights under unstable ambient conditions. In this program, the scale calculates the weight as the average value from a defined number of individual weighing operations. These weighing operations are also known as "subweighing operations" or "subweighs."

Available Features

- You can set the number of subweighing operations performed in the operating menu, or before starting the Averaging application (press [FUNCTION 2])
- The number of subweighs remaining to be performed is indicated in the display during weighing
- The measured result displayed is the arithmetic mean shown in the selected weight unit; a triangle under the plus or minus sign indicates that this is a calculated value
- Press [FUNCTION 1] for at least 2 sec. to display the pre-set number of subweighing operations
- Press [FUNCTION 2] to toggle the display between the measured result and the weight
- The measurement result is automatically output via the data interface port after initialization, if the menu code for "Printout with data ID codes" is set

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Preparation

Configure the Averaging application in the operating menu:

- If the scale is on, turn it off: Press [ON/OFF]
 - Turn on the scale: Press [ON/OFF]; while all segments are lit, press [TARE]
 - Select the “Averaging” menu item:
Press [TARE], [PRINT], and [PRINT]; press [TARE] repeatedly, if necessary
2 1 i2 Averaging
 - Confirm selection of Averaging: Press [PRINT]
for at least 2 sec.
 - > 0 indicates that this is the current menu code setting
 - Set the next parameters: Press [PRINT]
 - Select and confirm:
 - Number of subweighs for averaging:
 - 3 3 1 5 subweighs
 - 3 3 2 10 subweighs
 - 3 3 3 20 subweighs
 - 3 3 4 50 subweighs
 - 3 3 5 100 subweighs
- see also the “Scale Operating Menu (Overview)” in the chapter entitled “Configuring the Scale”
- Save settings and exit the Setup menu:
Press [TARE] for at least 2 sec.

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- Waage ausschalten (Taste [ON/OFF])
- Waage nullstellen (Taste [ZERO])
- Waage tarieren (Taste [TARE])
- Drucken (Taste [PRINT])

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- Start measuring ([FUNCTION 1] key)
- Clear display of results ([CAL/CF] key; press for at least 2 sec.)
- Change number of measurements after clearing the display ([FUNCTION 2] key)
- Display weight ([FUNCTION 2] key)
- Calibration/adjustment ([CAL/CF] key)

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Example

Determine the weight of a sample in extremely unstable ambient conditions. Select the number of subweighs while in the weighing mode.

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Menu: Application program: Averaging (2 / 12)

Step	Key (or instruction)	Display/Output
1. Turn on the scale	[ON/OFF]	
2. If necessary, zero the scale	[ZERO]	0.000 kg [#]
3. Set the number of subweighs for averaging (here: 5 subweighs)	[FUNCTION 2]	rEF 20 (briefly)
	[FUNCTION 2]	rEF 50 (briefly)
	[FUNCTION 2]	rEF 100 (briefly)
	[FUNCTION 2]	rEF 1 (briefly)
	[FUNCTION 2]	rEF 2 (briefly)
	[FUNCTION 2]	rEF 5 (briefly)
		0.000 kg ⁵
4. Place sample on the scale (weight readout fluctuates; here: around 275 g)		+ 8888 ⁵
5. Start measurement; if the print format is set to include data ID codes, the following is printed	[FUNCTION 1]	‡8888 ⁵ ... †
		‡0.275 kg ⁵
		Res + 0.275 kg
6. Unload the scale		‡0.275 kg ⁵
7. Repeat the procedure starting from Step 4.		

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Data Output Functions

There are 3 options for data output:

- Output to the display and control unit
- Output to a printer (generate a printout)
- Output to a peripheral device (e.g., computer) via the interface port

Printing a Data Record

Purpose

You can generate a printout of weights as well as other measured values and identifiers for documentation purposes. You can format the printout to meet individual requirements.

Available Features

The printout can be configured to include 2 user-defined header lines of 14 characters each. These text lines are entered either at the SCALTEC factory or on location with a special program that uses the BPI interface command sequence.

Print an individual weight, or net, gross and tare weights.

Line format: You can configure a data ID code of up to 6 characters for each of the values printed; this data ID code is printed at the beginning of the line

Printouts are generated automatically or by pressing the [PRINT] key, dependent on or independent of scale stability

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You can have the following values output automatically when using the application programs if menu code 7 1 2 or 7 1 3 is configured (printout with data ID codes):

- Second tare memory: Last net value
- Counting: Reference weight for one piece (average piece weight)
- Weighing in percent: Reference weight for the percentage selected
- Totalizing: Current weight, total weight (net)
- Averaging: Result of measurement

Factory Settings

Header:

The default header lines do not contain any text

Print manual/automatic:

Individual printout or automatic printout dependent on scale stability:

Manual after stability (5 1 2)

Line format:

Up to 6 characters at the beginning of each line to identify the weight or calculated value:

Print net, tare, gross value with data ID (7 1 3)

- See the chapter entitled "Configuring the Scale" for instructions on setting parameters

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Examples				
You can configure the printout to include 2 user-defined header lines. The software required for configuring these lines via your PC is available from your local SCALTEC office or dealer.	EISENMUELLER		User-defined	
	KASSEL		User-defined	
Printout without Data ID Codes:				
The value currently displayed is printed (weight or calculated value with unit)	+	1530.0 g	Weight in grams	
	+	58.562 ozt	Weight in Troy ounces	
	+	253 pcs	Piece count	
	+	88.2 %	Percentage	
	+	105.8 o	Calculated value	
Printout with Data ID Codes:				
The current value displayed can be printed with a data ID code of up to 6 characters at the beginning of the line. You can use this data ID code, e.g., to designate a weight readout as a net weight (N) or a calculated value as a piece count (QNT).	N	+	1.530 kg	Current net weight
	N1	+	1.530 kg	Current net weight
	T	+	0.234 kg	Value in tare memory
	T1	+	0.102 kg	Value in 2nd tare memory
	G	+	1.553 kg	Current gross weight
	G#	+	1.630 kg	Calculated gross weight
	Qnt	+	253 pcs	Calculated quantity
	Prc	+	88.2 %	Calculated percentage
	Sum	+	1.279 kg	Total value (net)
	Res	+	1.530 kg	Calculated result

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Print Application Parameters:

You can generate a print-out of one or more of the values configured for initialization of an application as soon as you initialize the scale.

wRef + **0.014 kg**
wxx% + **1.200 kg**

Counting:
Average piece weight
Weighing in percent:
reference weight for the selected percentage

Auto Print:

You can have the weight readout printed automatically. The display update interval depends on the operating status of the scale and on the scale model.

N + **1.530 kg**
stat
stat L
stat H

Net weight
Display blank
Display underload
Display overload

For details on the data interface (such as the data output or input formats, pin assignment, etc.), contact SCALTEC.

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Error codes are shown in the main display for approx. 2 seconds, after which the program automatically returns to the weighing mode.

Display	Cause	Solution
No segments appear on the display	No AC power is available	Check the AC power supply
	The AC adapter is not plugged in	Plug in the AC adapter
	Automatic shutoff is configured	Turn on the scale
	Rechargeable battery has run down	Recharge battery (see battery instructions)
H	The load exceeds the scale capacity	Unload the scale
L	The load plate is not in place	Place the load plate on the scale
	Something is touching the load plate	Move the object that is touching the load plate
E 01	Data output not compatible with input format	Change the configuration in the Setup menu
E 02	Calibration parameter not met; e.g.:	Calibrate only when zero is displayed
	– scale not zeroed	Press [ZERO] to zero the scale
	– scale is loaded	Unload the scale

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Display	Cause	Solution
E 08	The scale was zeroed outside the zero range	Only zero the scale when in the zero range; i.e.: (2% of the maximum capacity)
E 09	Taring is not possible when the gross weight is \leq zero	Press [ZERO] to zero the scale
E 10	The [TARE] key is blocked when there is data in the second tare memory (net-total) – only 1 tare function can be used at a time	Press [CAL/CF] to clear the tare memory and release the tare key
E 11	Tare memory not accessible	Press [ZERO]
E 22	Weight is too light or there is no sample on the scale	Increase the reference quantity
E 30	Interface port for printer output is blocked	Perform "Reset menu" or Contact your local SCALTEC Service Center
The maximum weighing capacity is less than is indicated in "Specifications"	The scale was turned on without the load plate in place	Press the [ON/OFF] key to turn the scale off and back on again

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Display	Cause	Solution
The weight readout changes constantly	Unstable ambient conditions	Set up the scale in another area
	Too much vibration, or the scale is exposed to a draft	Change setup configuration to adapt the scale to the ambient conditions
The weight readout is obviously wrong	A foreign object is caught between the load plate and the scale frame	Remove the foreign object
	The scale has not been calibrated/adjusted	Calibrate/adjust the scale
	The scale was not tared before weighing	Tare before weighing

If any other errors occur, contact your local SCALTEC Service Center!

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Service

Regular servicing by a SCALTEC technician will extend the service life of your scale and ensure its continued weighing accuracy.

Repairs

Repair work must be performed by trained service technicians. Any attempt by untrained persons to perform repairs may lead to hazards for the user.

Cleaning

- ⚠ Disconnect the scale from the AC adapter and unplug any data cables that are connected to the scale
- ⚠ Make sure that no liquids enter the scale housing
- ⚠ Do not use any aggressive cleaning agents (solvents or similar agents)
- Unplug the AC adapter from the wall outlet (mains supply)
- If you have a data cable connected to the interface, unplug it from the scale
- Clean the scale using a piece of cloth which has been wet with a mild detergent (soap)
- After cleaning, wipe down the scale with a soft, dry cloth

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Safety Inspection

If there is any indication that safe operation of the scale with the AC adapter is no longer warranted:

- Turn off the power and disconnect the equipment from AC power immediately
- > Lock the equipment in a secure place to ensure that it cannot be used for the time being

Safe operation of the scale with the AC adapter is no longer ensured when:

- there is visible damage to the AC adapter
- the AC adapter no longer functions properly
- the AC adapter has been stored for a relatively long period under unfavorable conditions

In this case, notify your nearest SCALTEC Service Center. Maintenance and repair work may only be performed by service technicians who are authorized by SCALTEC and who

- have access to the required maintenance manuals
- have attended the relevant service training courses

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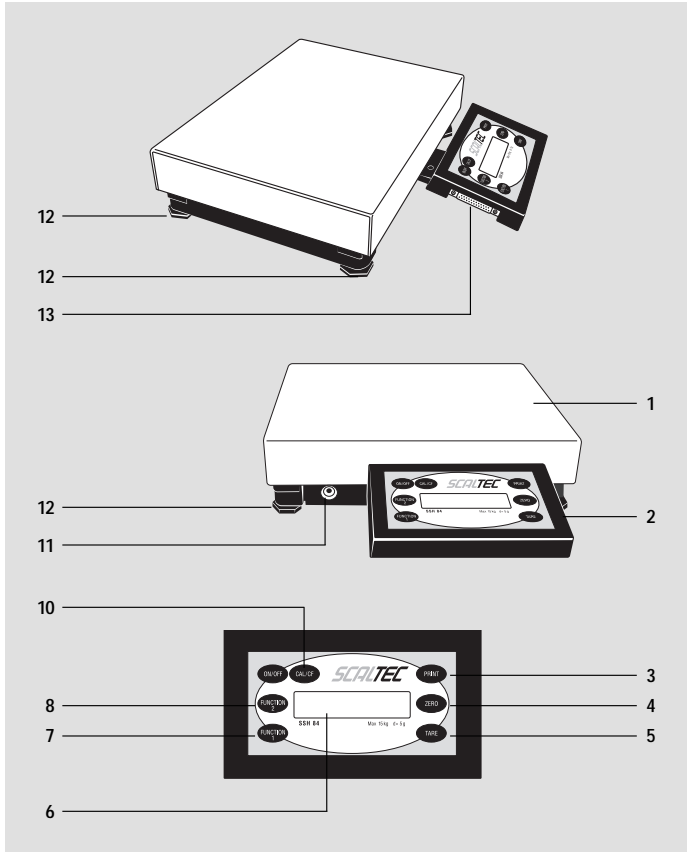
Instructions for Recycling the Packaging

To ensure safe shipment, your scale has been packaged using environmentally friendly materials. After successful installation of the scale, you should recycle this packaging.

For information on recycling options, including recycling of old weighing equipment, contact your municipal waste disposal center or local recycling depot.

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¹⁾ to delete: press the press the "CAL/CF" key for more than 2 sec.

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Description of the Keys

[ON/OFF] key: On/off switch

Switches the display on/off. The scale remains in the standby mode.

[ZERO] key: Zero

Sets the readout to zero. The scale can only be zeroed if the load is $\pm 2\%$ of the maximum capacity.

[TARE] key: Tare

For taring the weight of containers so that the scale readout shows the net weight of samples.

[CAL/CF] key: Calibrate/adjust or Clear function

Press briefly (under 2 sec.) to:

- Start the calibration function

Press for at least 2 seconds to:

- Clear memory and delete application data
- Interrupt calibration/adjustment routines

[FUNCTION 1] key: Start Application Program

The procedure that follows after this key is pressed depends on the application program selected; for a complete description, see the corresponding section in the chapter entitled "Operating the Scale".

[FUNCTION 2] key: Toggle

For toggling the display between weighed and calculated values (counting result, percentage, calculated result)

For changing the reference quantity/percentage or numbers of subweighs for averaging

[PRINT] key: Data Output

Press this key to output the displayed value via the built-in data interface port to a SCALTEC "Data Printer" or a computer.

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Standard models

Model		SSH84	SSH83	SSH82	SSH81
Weighing capacity	kg	15	35	60	150
Readability	g	5	10	20	50
Maximum load	kg	30	70	120	300
Tare range (subtractive)	kg	15	35	60	150
Repeatability	≤+g	5	10	20	50
Linearity	≤+g	5	10	20	50
External calibration weight (of at least accuracy class...)	kg lb	5 (M1) 10 (M1)	10 (M1) 20 (M1)	20 (M1) 50 (M1)	50 (M1) 100 (M1)
Net weight, approx.	kg	5	8	24	24
Dust and water protection rating for the weighing platform according to EN 60529		IP54			
Dust protection rating for the display and control unit according to EN 60529		IP40			
AC power source/power requirements		AC adapter, 230 or 115 V, +15% ... -20%			
Frequency		48 – 60 Hz			
Power consumption		maximum 16 VA; average 8 VA			
Operating temperature range		-10 ... +40 °C (263 ... 313 K, 32 °F ... 104 °F)			
Adaptation to ambient conditions		By selection of 1 of 4 optimized filter levels			
Display update (depends on the filter level selected)		0.1 – 0.4			
Hours of operation with fully charged SBP01 external battery pack, approx.		48 h			
Selectable weight units		Grams, kilograms, carats, pounds, ounces, Troy ounces, Hong Kong taels, Singapore taels, Taiwanese taels, grains, pennyweights, milligrams, parts per pound, Chinese taels, mommes, Austrian carats, tola, baht, mesghal and tons			
Built-in interface		RS-232			
Format:		7-bit ASCII, 1 start bit, 1 or 2 stop bits			
Parity:		mark, odd, even or space			
Transmission rates:		150 to 9,600 baud			
Handshake mode:		Software or hardware			

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
Specifications

Model		SSH94	SSH93	SSH92	SSH91
Weighing capacity	kg	15	35	60	150
Readability	g	0.5	1	2	5
Maximum load	kg	30	70	120	300
Tare range (subtractive)	kg	15	35	60	150
Repeatability	≤±g	0.5	1	2	5
Linearity	≤±g	1	2	4	10
External calibration weight (of at least accuracy class...)	kg lb	5 (F2) 10 (F2)	10 (F2) 20 (F2)	20 (F2) 50 (F2)	50 (F2) 100 (F2)
Net weight, approx.	kg	5	8	24	24
Dust and water protection rating for the housing weighing platform according to EN 60529	IP54				
Dust protection rating for the housing weighing platform according to EN 60529	IP40				
AC power source/power requirements	AC adapter, 230 or 115 V, +15% ... – 20%				
Frequency	48 – 60 Hz				
Power consumption	maximum 16 VA; average 8 VA				
Operating temperature range	–10 ... +40 °C (263 ... 313 K, 32 °F ... 104 °F)				
Adaptation to ambient conditions	By selection of 1 of 4 optimized filter levels				
Display update (depends on the filter level selected)	0,1 – 0,4				
Hours of operation with fully charged SBP01 external battery pack, approx.	48 h				
Selectable weight units	Grams, kilograms, carats, pounds, ounces, Troy ounces, Hong Kong taels, Singapore taels, Taiwanese taels, grains, pennyweights, milligrams, parts per pound, Chinese taels, momes, Austrian carats, tola, baht, mesghal and tons				
Built-in interface	RS-232				
Format:	7-bit ASCII, 1 start bit, 1 or 2 stop bits				
Parity:	mark, odd, even or space				
Transmission rates:	150 to 9,600 baud				
Handshake mode:	Software or hardware				

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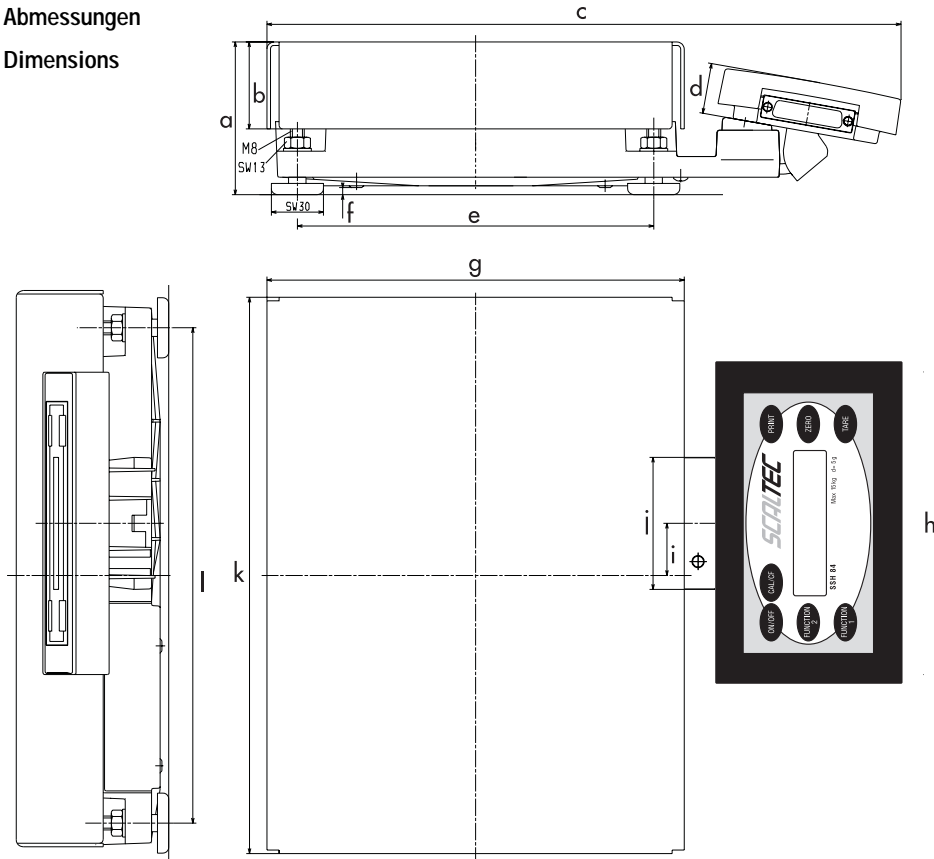Data Printer

SPR01

with date/time, statistical data
evaluation and transaction counter
functions (can only be operated
with an AC adapter)

AC adapter for the Data Printer for use in

Europe

629801

USA

629802

Australia

629804

Paper (5 rolls)

SPR01-1

External rechargeable battery pack

SBP01

has a battery-level indicator (LED); can be recharged using the
AC adapter (time it takes to charge the discharged battery pack:
15 hours); see "Specifications" for hours of operation

Calibration weights

**Information
on request**

for all SSH scales; extensive assortment, optionally
available with officially recognized DKD certificates (DKD
is the German Calibration Service)

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The CE marking affixed to the equipment indicates that the equipment meets the requirements of the following Directive(s) issued by the Council of the European Union:

Council Directive 89/336/EU "Electromagnetic compatibility (EMC)"

Important Note:

The operator shall be responsible for any modifications to SCALTEC equipment and for any connections of cables or equipment not supplied by SCALTEC and must check and, if necessary, correct these modifications and connections. On request, SCALTEC will provide information on the minimum operating specifications (in accordance with the Standards listed above for defined immunity to interference).

73/23/EU "Electrical equipment designed for use within certain voltage limits"

If you use electrical equipment in installations and under ambient conditions requiring higher safety standards, you must comply with the provisions as specified in the applicable regulations for installations in your country.

For information on the current legal requirements of your country, please contact your local SCALTEC customer service office.

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Index	
Accessories	Page 70
Application programs	Page 29
Auto Print	Page 55
Auto zero	Page 20
Averaging	Page 51
Basic weighing function	Page 24
Calculation by a factor	Page 43
Calibration/adjustment	Page 27
Care and maintenance	Page 62
CE Marking	Page 72
Cleaning the scale	Page 62
Configuring the scale	Page 14
Connecting AC power	Page 8
Contents	Page 3
Counting	Page 29
Data output functions	Page 55
Description of the keys	Page 65
Dimensions	Page 68
Display	Page 13
Equipment supplied	Page 6
Error codes	Page 59
Factory settings	Page 18
Filter adaptation	Page 20
General view of the scale	Page 64
Getting started	Page 5
Gross weight: toggling	Page 49
Installation instructions	Page 7
Keys: description of	Page 65
Leveling the scale	Page 11
Menu (overview)	Page 20
Net weight: toggling	Page 49
Net-total (2nd tare memory)	Page 36
Operating design	Page 12
Operating menu (overview)	Page 20
Operating the scale	Page 24
Options (Accessories)	Page 70
Printing	Page 55
Printing: manual/automatic	Page 56
Recycling the packaging	Page 63
Repairs	Page 62
Safety inspection	Page 63
Safety precautions	Page 4
Setting parameters in the menu	Page 17
Specifications	Page 66
Stability range	Page 20
Tare memory	Page 41
Taring	Page 24, 26
Toggle:	
between gross/net weights	Page 49
between weight units	Page 32
Totalizing	Page 46
Unit indicator	Page 13
Unpacking the scale	Page 6
Warmup time	Page 10
Warranty	Page 5
Weighing in percent	Page 40
Weight unit: toggling	Page 32
Weight units	Page 34
Zeroing the scale	Page 24, 26

**SCALTEC
Instruments GmbH**

Rudolf-Diesel-Strasse 1
37308 Heiligenstadt, Germany
Phone (3606) 6701-0
Telefax (3606) 6701-60
E-Mail: Scaltecba1@ad.com

Copyright by Scaltec Instruments GmbH, Heiligenstadt, Germany.
All rights reserved.

No part of this publication may be reprinted or translated in
any form or by any means without the prior written permission
of Scaltec Instruments GmbH.

The status of the information, specifications and illustrations
in this manual is indicated by the date given below.

Scaltec Instruments GmbH reserves the right to make changes to the technology,
features, specifications, and design of the equipment without notice.

Status: July 1998, Scaltec Instruments GmbH, Heiligenstadt, Germany.

98648-007-93

Printed in Germany
on paper that has been bleached without any use
of chlorine · JW/ZPD · W398-A52.SSH · M
Publication No. WXX6034-a98071

SCALTEC